

Lab Leg Panel


Product Numbers

- LULPxyyy


Required Tools

- 5mm Allen Wrench


Hardware Pack: LULPHP (p/n 09061701-01)

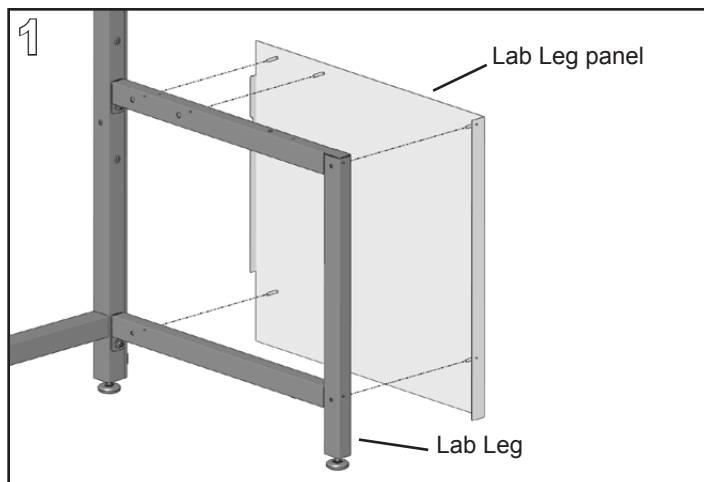

Align the PEM studs in the Lab Leg Panel with the holes in the side of the Lab Leg.

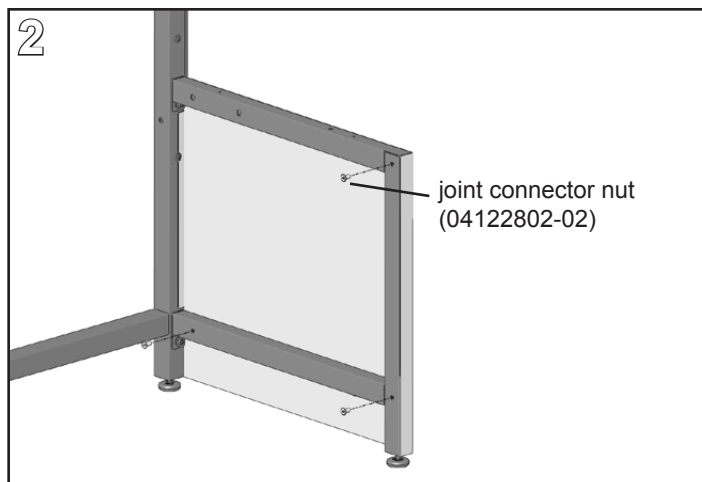

Loosely fasten the panel to the leg with a joint connector nut through the bottom leg holes and the top front hole in the leg.

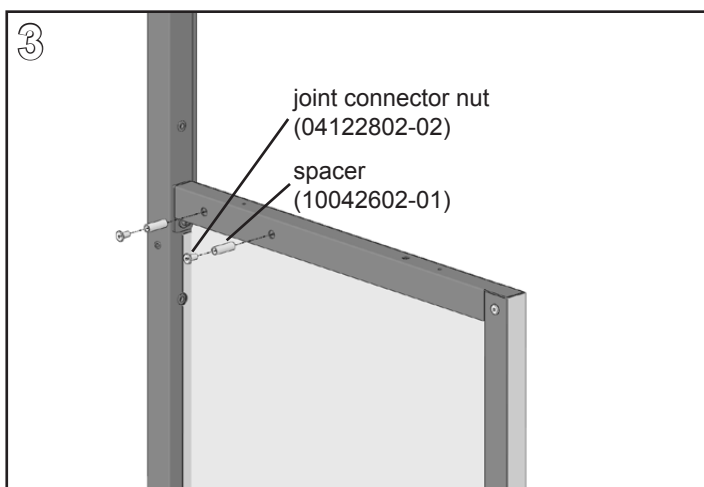

Insert a spacer into each of the holes in the top channel of the leg (24" deep has one hole, 30" and 36" depths have two holes). Fasten with a joint connector nut through each spacer/leg/bracket assembly. Tighten all nuts securely.

SYMBIOTE® Technical Furniture

www.symbiote.com 300 N. Centennial Street Zeeland, MI 49464 phone 616.772.1790 fax 616.772.1797

p/n 09021216-01

sheet 1 of 1